


# MOBILE LAB UNITS

4111, 4121, 4222, 4332, 4342, & 4352 SERIES (OAK OR MAPLE)  
ASSEMBLY INSTRUCTIONS

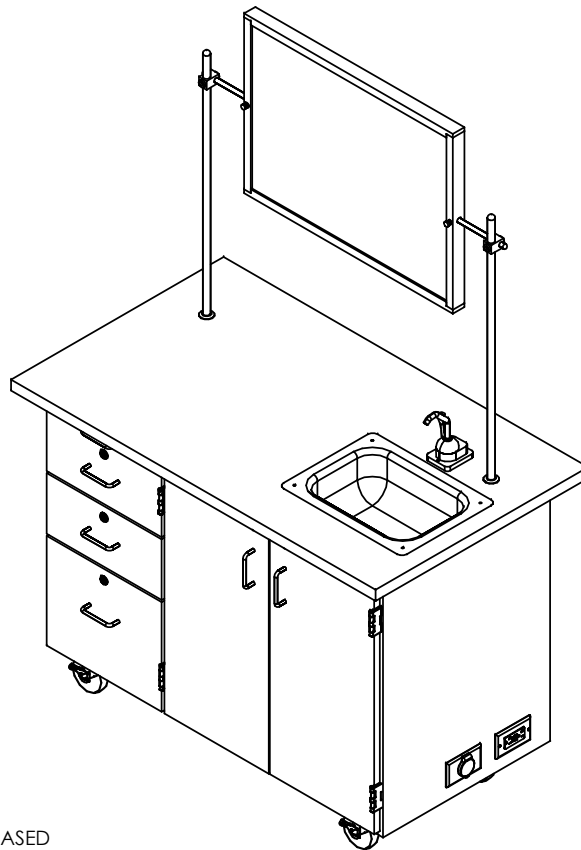

NOTE:

1. MODEL VIEWS MAY NOT REPRESENT EXACT MODEL PURCHASED



[StoreMoreStore.com](http://StoreMoreStore.com) | 855-786-7667

## TOOLS REQUIRED


½" WRENCH
ADJUSTABLE PLIERS
SCREW DRIVER

## ASSEMBLY COMPONENTS

ITEMS INCLUDED	PART #	PART DESCRIPTION	QTY
CASTERS	100081	CASTER,4IN SWIVEL TOP PLATE, W/BRAKE	4
CASTER HARDWARE KIT	206509	CASTER,ASSY,HDWRE,16-SCREWS, 16-NUTS	1
**EXTENSION CORD**	100036	ELECTRICAL EXTENSION CORD	1
**HAND PUMP**	100013	PUMP,ROCKET HAND RP800	1
**TUBING**	100015	TUBING,⅜ FOR HAND PUMP	2.5
**CROSS BAR**	100007	ROD,CROSSBAR ½ DIA.X 40L	1
**UPRIGHTS**	100004	ROD,UPRIGHT ¾ DIA.X 40L	2
**CLAMPS**	100003	ROD CLAMP ½X¾ ASSY	2
**WATER JUGS**	100118	WATER JUG,2 1/2 GAL RECTANGULAR	2
**GOOSE NECK FIXTURE**	222830	FIXTURE,L081-55WSA GSNCK W/ FLANGE	1

### NOTE:

1. COMPONENTS VARY WITH STYLE OF MOBILE CABINET PURCHASED.
2. \*\*XX\*\* INDICATES ITEMS THAT MAY NOT COME WITH UNIT (DEPENDING ON MODEL PURCHASED).

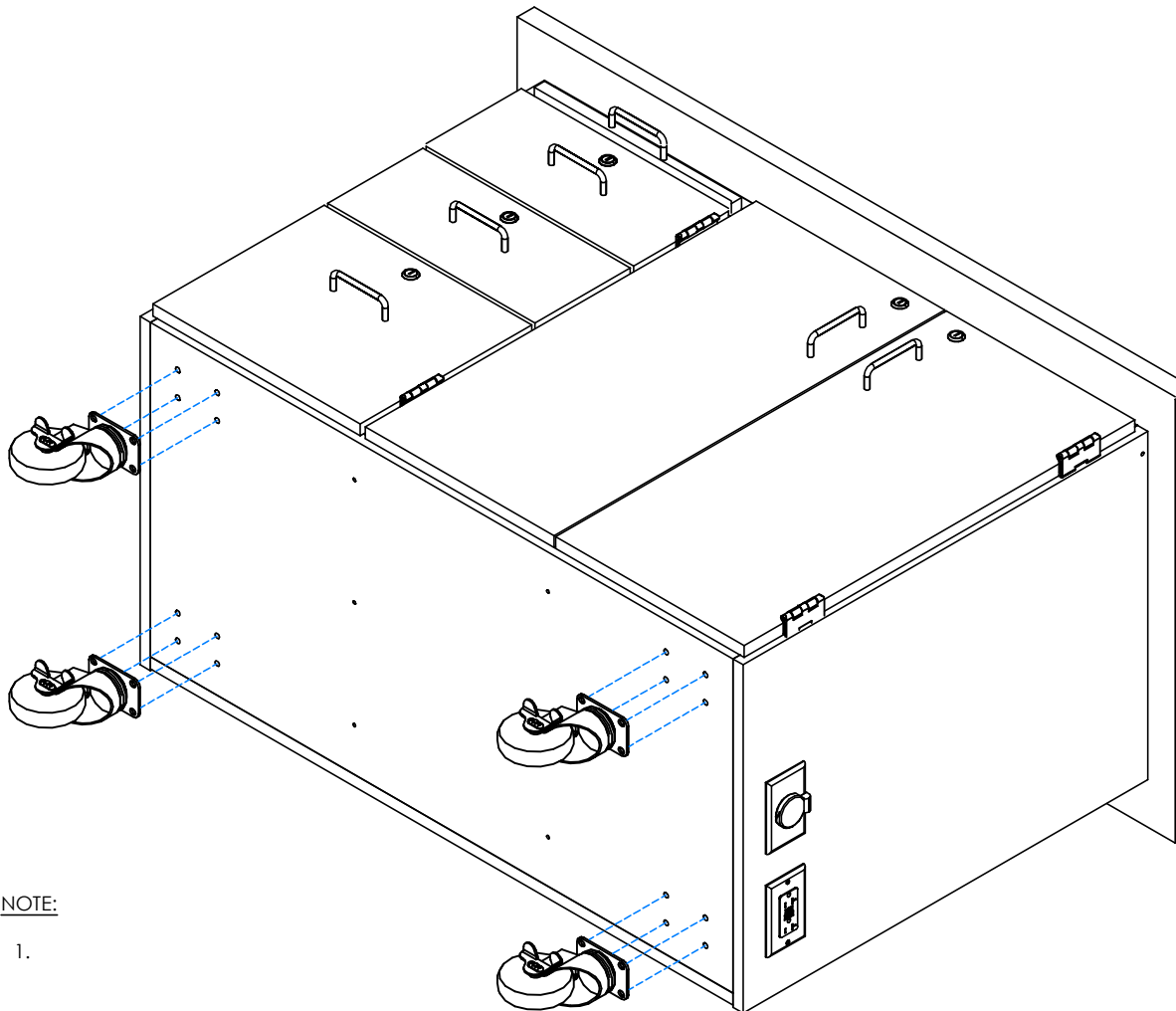


NOTE:

1.

STEP 1:


- LAY CABINET ON ITS BACK.
- ATTACH CASTERS FROM INSIDE OF CABINET USING HARDWARE PROVIDED.
- IF YOUR CABINET DOES NOT HAVE A HAND PUMP, WATER FIXTURE, OR UPRIGHTS AND A CROSSBAR, SKIP TO STEP 6.

**\*\*NOTE\*\***  
BE SURE CABINET IS ON  
A CLEAN SURFACE TO  
PREVENT DAMAGE

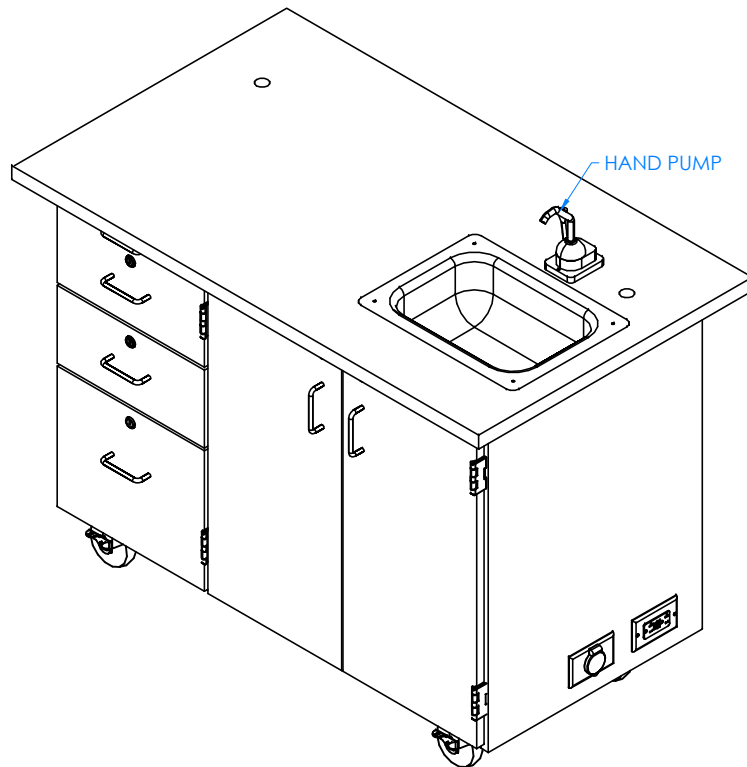

4332 UNIT SHOWN WITH HAND PUMP

GOOSE NECK FIXTURE

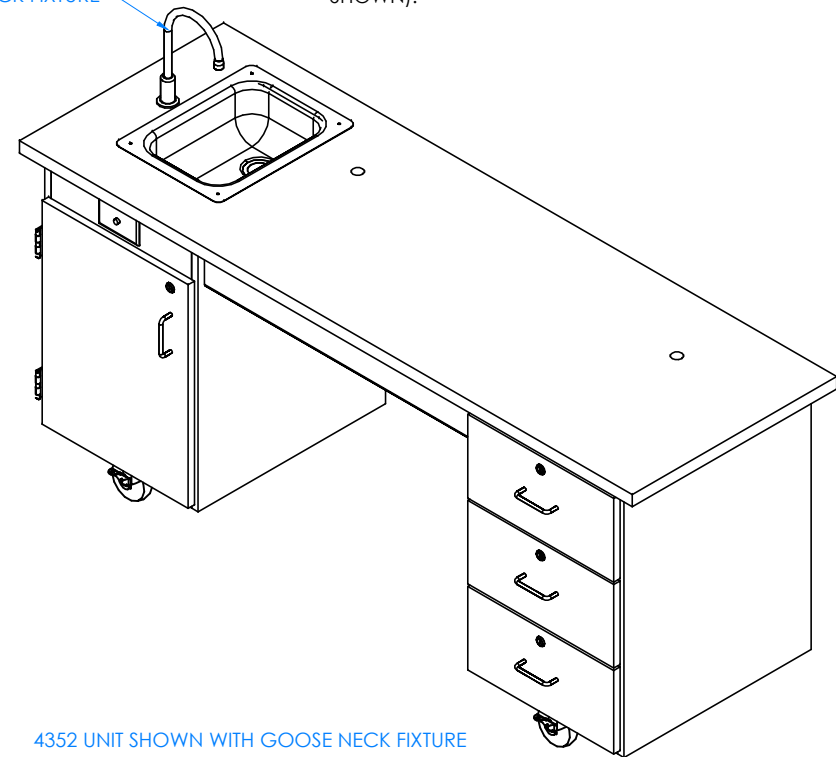


4352 UNIT SHOWN WITH GOOSE NECK FIXTURE

STEP 2:

- ATTACH HAND PUMP OR WATER FIXTURE TO PILOT HOLES LOCATED ON TOP WITH HARDWARE PROVIDED.
- ATTACH ONE END OF THE  $\frac{3}{8}$ " TUBING TO THE PUMP OR WATER FIXTURE AND INSERT OTHER END INTO FRESH WATER CONTAINER (NOT SHOWN).
- REMOVE SINK STOPPER FROM BOTTOM OF PLASTIC SINK DRAIN HOSE AND PUT INTO THE SINK DRAIN.
- INSERT END OF THE HOSE INTO WASTE WATER CONTAINER (NOT SHOWN).

NOTE:

1.

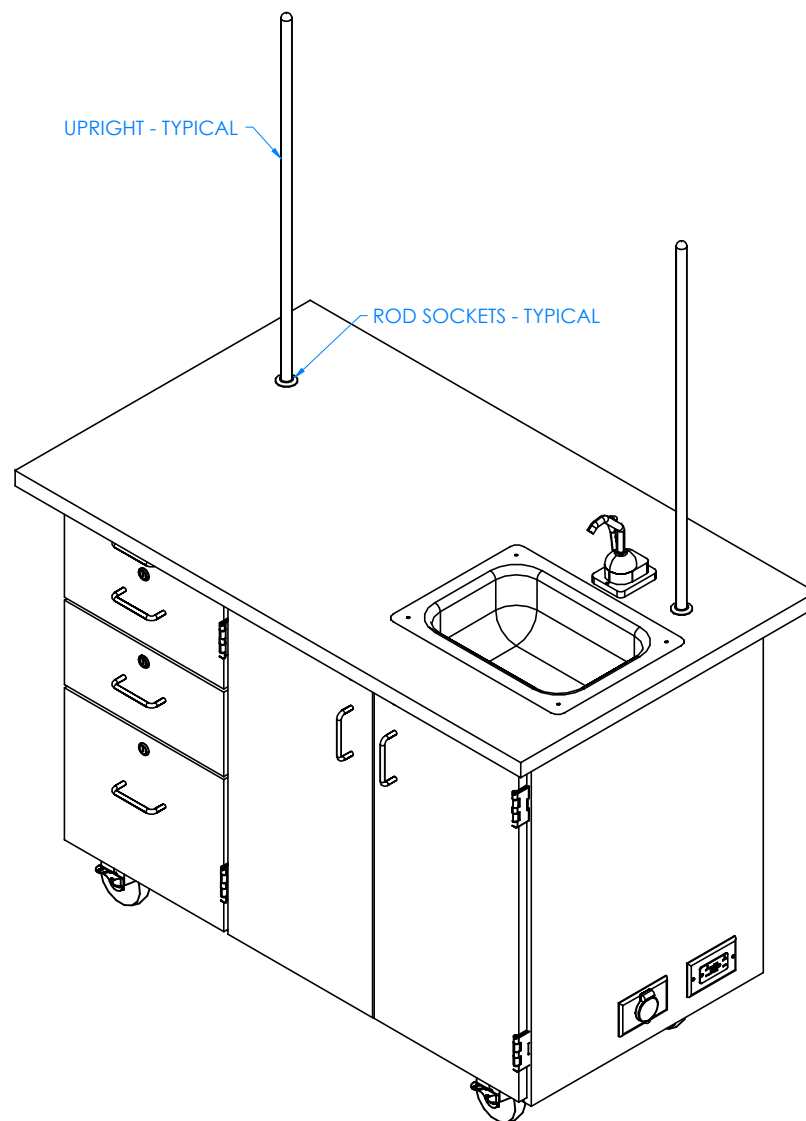


STEP 3:

- INSERT TAPERED END ON UPRIGHTS INTO ROD SOCKETS ON THE TOP.

NOTE:

1.

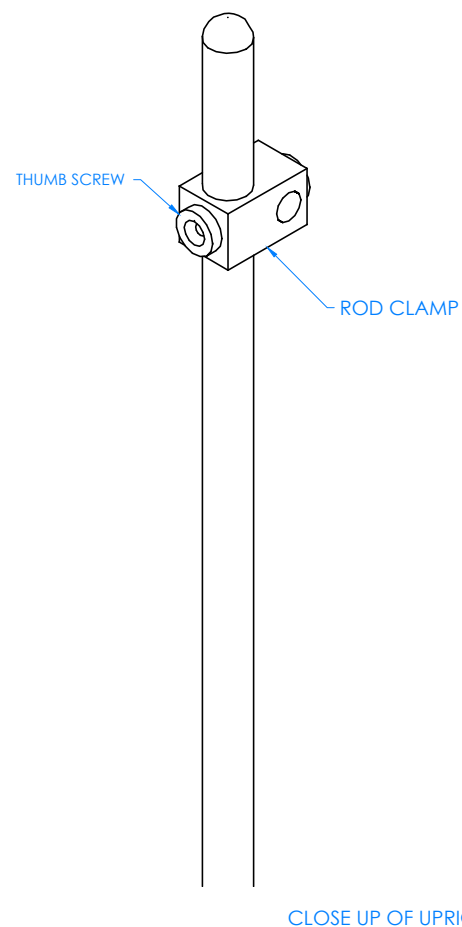


STEP 4:

- SLIDE ROD CLAMPS OVER UPRIGHT AND TIGHTEN IN DESIRED LOCATION USING THUMB SCREWS PROVIDED ON CLAMPS.
- REPEAT ON OTHER UPRIGHT.

NOTE:

1.

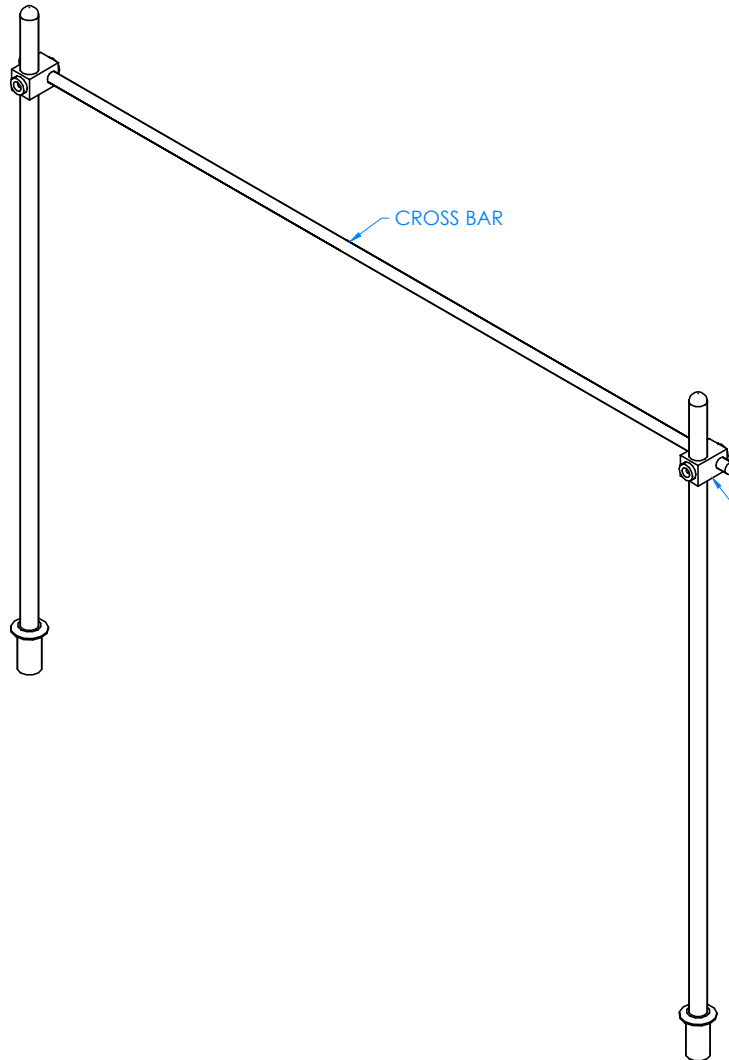


STEP 5:

- SLIDE CROSS BAR THROUGH ONE CLAMP.
- POSITION OTHER SIDE THROUGH THE OTHER CLAMP.
- IF YOU HAVE A MIRROR/MARKER BOARD, SEE MIRROR/MARKER BOARD ASSEMBLY INSTRUCTIONS FOR PROPER INSTALLMENT.

NOTE:

1.

CLOSE UP OF UPRIGHT AND  
CROSSBAR ASSEMBLY

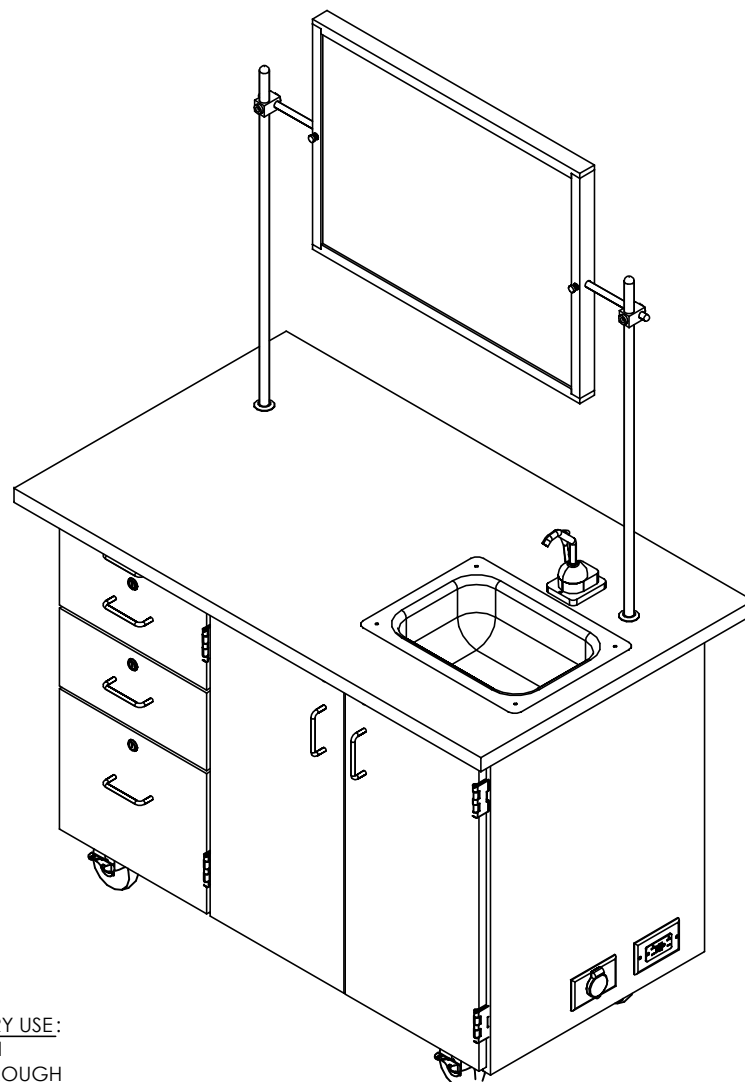

STEP 6:

- SET CABINET IN DESIRED LOCATION.

NOTE:

1. RECOMMENDED CLEANING INSTRUCTIONS AFTER EVERY USE:  
EMPTY CLEAN AND WASTE JUGS. RINSE WITH MILD ANTI BACTERIAL SOAP. RUN MILD ANTI BACTERIAL SOAP THROUGH WATER LINES AND HAND PUMP TO INSURE NO BACTERIAL BUILD UP.

REV.

COMPLETE UNIT SHOWN